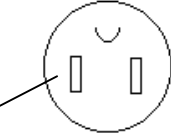
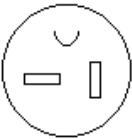
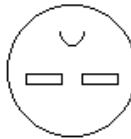



## Plugs For HA And HV Series Autoclaves (View With Pins of Plugs Facing You)

	HVE-25/HV-25	HVE-50/HV-50 HVP-50/HA-300	HV-85	HV-110
Connect white neutral wire to this pin.				
Plug type:	125V, 15A NEMA 5-15P	125V, 20A NEMA 5-20P	240V, 15A NEMA 6-15P	240V, 20A NEMA 6-20P
Heater:	1.5 KW(1500VA)	2.0 KW(2000VA)	3.0 KW(3000VA)	4.0 KW(4000VA)
Phase:	SINGLE	SINGLE	SINGLE	SINGLE

**NOTE:** You can operate HV-85 and HV-110 autoclaves from a 208V, 220V, or 240V power source.

If the outlet can support the amperage required by the heater, but the plug does not fit the outlet, have an electrician change the outlet to fit the plug or, the plug to fit the outlet. For example, a 240V, 30A outlet can support an HV-85 or an HV-110 and the autoclave plug can be replaced by a 240V, 30A plug.