


Front Panel Components

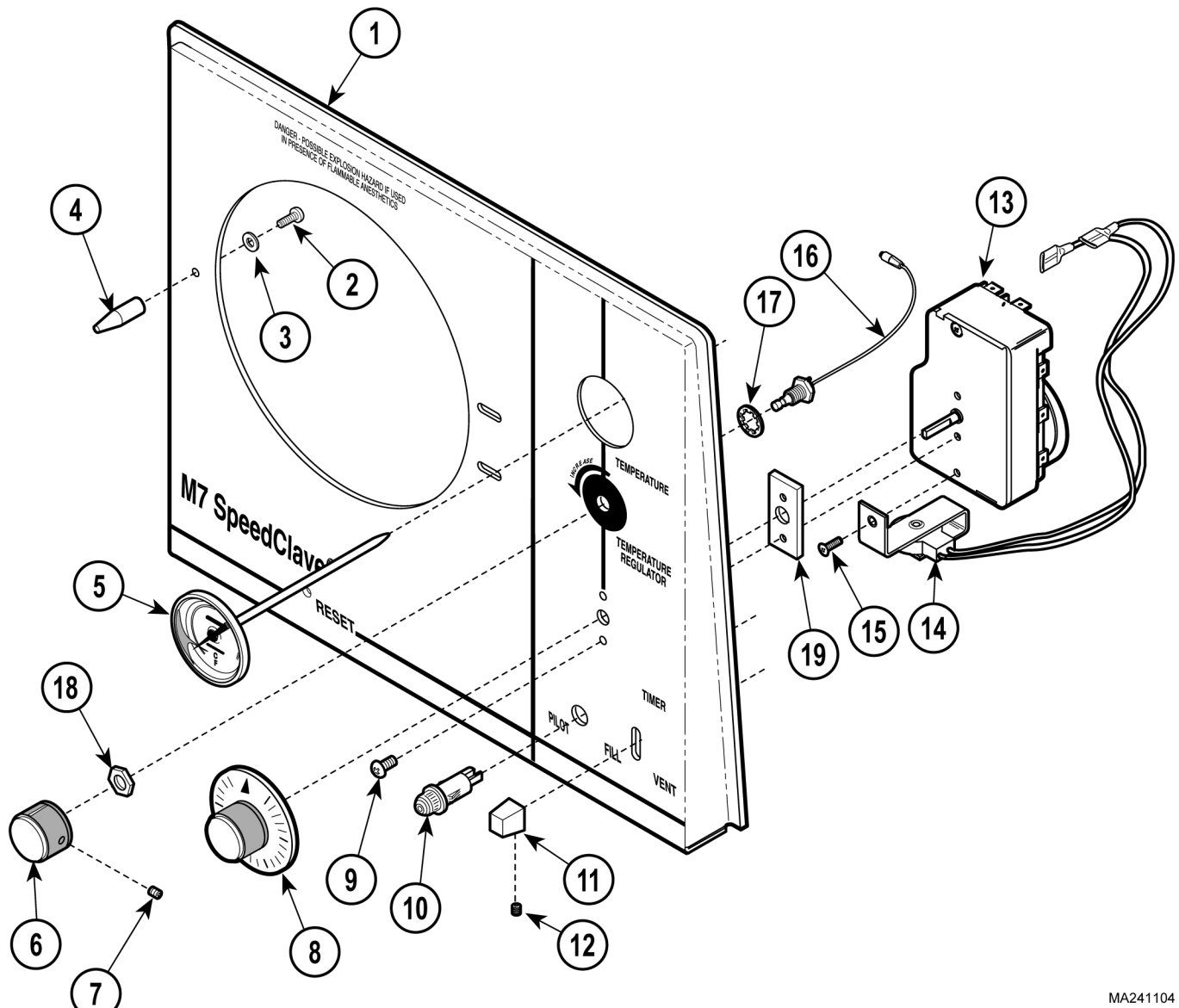

MA241104

**Used on Units with Serial Numbers RB18475 thru RB21249,
CR1725 thru CR1904, CS8319 and CS11602, FM1546 thru FM1875, and FN1310
thru FN1499**

Item	Part No.	Description	Qty.	Item	Part No.	Description	Qty.
1	MZY100067 MZY100068	Ritter Panel (M7-001) Midmark Panel (M7-002, M7-003, M7-004)	1	11	MZY100073	Fill / Vent Knob Assembly (Includes Item 12)	1
	MZY100069	Special Panel (M7-005, M7-006)	1	12	• MZY100058	• Setscrew	1
2	MZY100075	Screw	1	13	MZY100059	Timer (M7-001, M7-003, M7-004, M7-005) [Incl. Items 14 and 15]	1
3	MZY100071	Lockwasher	1		MZY100060	Timer (M7-002, M7-006) [Incl. Items 14 and 15]	1
4	MZY100076	Alignment Pin	1		MZY100077	Timer (Units with "DR" S/N Prefix)	1
5	MZY100052	Thermometer	1	14	• MZY100064	• Timer Buzzer (120 volt)	1
6	MZY100053	Control Knob (Includes Item 7)	1		• MZY100065	• Timer Buzzer (230 volt)	1
7	•	• Setscrew	1	15	• MZY100063	• Screw	1
8	MZY100054	Timer Knob	1	16	MZY100049	Flexible Shaft Assembly (Includes Item 17 and 18)	1
9	MZY100063	Screw	2	17	• MZY100050	• Lockwasher	1
10	MZY100055	Pilot Light (M7-001, M7-003, M7-004, M7-005)	1	18	• MZY100051	• Nut	1
	MZY100056	Pilot Light (M7-002, M7-006)	1	19	MZY100074	Spacer	1

Always Specify Model & Serial Number