


Tank Assembly

SECTION VI PARTS LIST

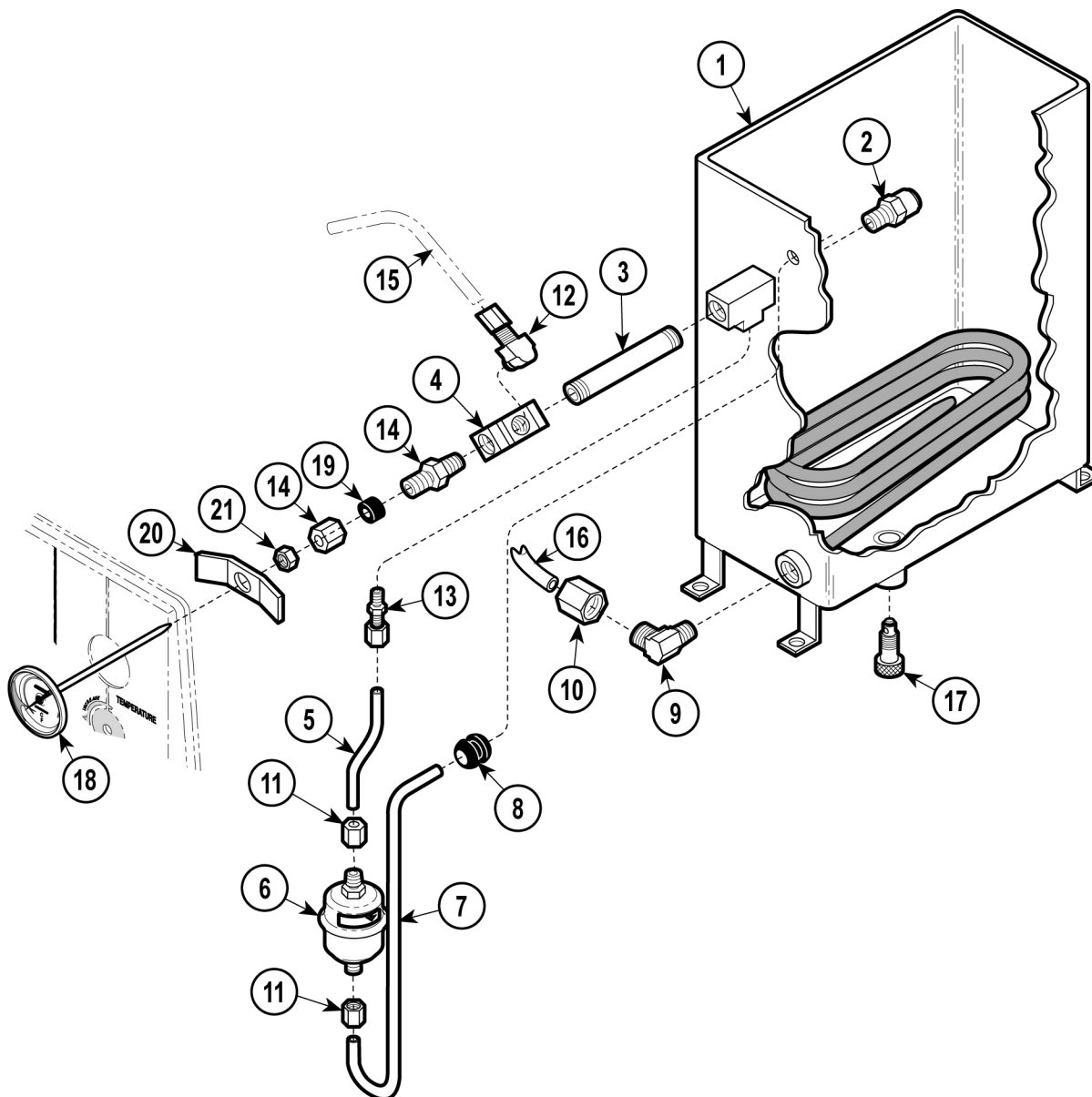

MA240808

Used on Units with Serial Numbers RB7987 thru RB17992, CR1158 thru CR1685, DR1000 thru DR1267, CP1062 thru CP1149, CS1976 thru CS7753 FM1000 thru FM1425 and FN1000 thru FN1229

Item	Part No.	Description	Qty.	Item	Part No.	Description	Qty.
	MZY100185	Tank Assembly (Incls. Items 1 thru 21)	1	12	MZY100175	• Male Connector	1
1	MZY100186	• Tank	1	13	MZY100135	• Connector Drilled	1
2	MZY100170	• Safety Valve (Apply Teflon Tape #014-0026-00)	1	14	MZY100179	• Compression Nut	1
3	MZY100171	• Pipe Nipple	1	15		Tube (Refer to "Sterilizer Main Components" Elsewhere)	Ref
4	MZY100172	• Pipe "T"	1	16		Tube (Refer to "Sterilizer Main Components" Item 23)	6-5.1
5	MZY100190	• Tube	1	17	MZY100182	• Drain (Includes O-ring)	1
6	MZY100174	• Bellows	1		MZY100183	• O-ring	1
7	MZY100191	• Tube	1	18	MZY100052	• Thermometer (Includes Item 26)	1
8	MZY100189	• Grommet	1	19	MZY100184	• Thermometer Sleeve	1
9	MZY100137	• Male Elbow	1	20	MZY100180	• Bracket	1
10	MZY100178	• Compression Nut	1	21	MZY100181	• Thermometer Nut	1
11	MZY100136	• Compression Nut	4				

Always Specify Model & Serial Number